



University of Melbourne iGEM 2014 Lab Procedure

Procedure	Name:	Cleanup		
	Version:	2		
	Description:	How to dispose of live cultures and clean the flasks Cleanup Video		
	Trigger:	Use this procedure when finished with live cultures		
Last updated	Name:	Robyn Esterbauer	Date:	22.06.14
You will need	Time:	10 minutes/overnight		
	PPE:	Gloves Safety glasses Lab coat		
	Equipment:	Fume hood		
	Materials:	Bleach		
Day 1				
Step 1	Take culture flask to fume hood			
Step 2	Transfer to plastic bottle			
Step 3	Add 100mL of bleach for every 1L of culture, leave overnight			
Day 2				
Step 4	Tip the solution left overnight down the drain in the fume hood			
Step 5	Rinse the bottle with water			
Step 6	Place the flask with the other items to be autoclaved with foil and autoclave tape on the top Note: for autoclave instructions see ‘			
Version history	Revised by Elizabeth Brookes on 16.07.14: <ul style="list-style-type: none"> • Use of hypo eliminated. 			