

## Blood Agar (D-Agar)

•

Glucose	1g
Bactopeptone	10g
Neopetone	5g
Yeast extract	1,25g
NaCl	5g
Tris	1,25g
H <sub>2</sub> O	ad1000ml

- autoclave
- cool down to 48 °C
- add 3 ml sheep blood per 100ml

Protocol generously provided by the lab  
Prof. Thorsten Mascher  
Großhadernerstr. 2-4  
82152 Planegg-Martinsried  
[www.syntheticmicrobe.bio.lmu.de](http://www.syntheticmicrobe.bio.lmu.de)